

Use of the Microscope

Please refer to the manuals in the microscope room or ask for assistance.

Powering on:

Make sure the Openlab program is closed

Turn on the microscope

Turn on the mercury light (if necessary)

Turn on the camera box

Cleanup:

If you used the Oil Immersion objective, clean it by pulling a Kimwipe over it

DO NOT WIPE OR APPLY STRONG PRESSURE AS THIS MAY DAMAGE THE LENS

Lower the stage

If you used the mercury lamp, record the counter number when you turn it off

THE LAMP MUST BE ON FOR AT LEAST 30MIN BEFORE IT CAN BE TURNED OFF