

Using Experiment Templates in SpectroFlo

Written by: Laura Johnston

Date: 07/06/2020

Revised: 09/21/2020


Data Management

- Data is pushed from the aurora computer to the server every 30 minutes.
- On the Aurora computer: Once a week, anything in the “My Experiments” window in SpectroFlo that is older than 30 days will be deleted from the software.
- The data uploaded on the server will be discarded after one year. The long term storage of data remains your own responsibility and you should develop your own data storage strategy. We recommend a cloud storage service, such as [UChicago box](#) or dropbox.


NOTE: Templates will not be deleted on a regular basis.

Create a template by saving an experiment as a template


1. If you would like to save your experiment template, go back to the acquisition tab, open your experiment and select Save As

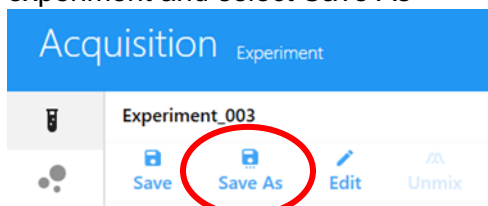

2. Name your template (including your name and date is recommended)
3. Your template can now be found in the template window

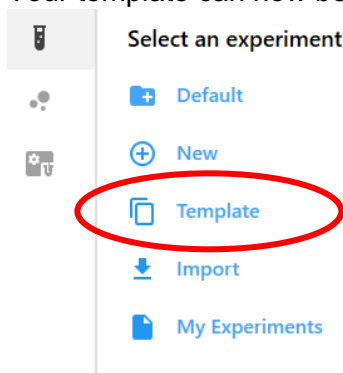


NOTE: After opening a new experiment from template, remember to rename the experiment with your name and date

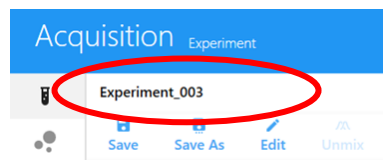

To export your saved template


1. Go to the library tab and select experiment template
2. Select your template and then export

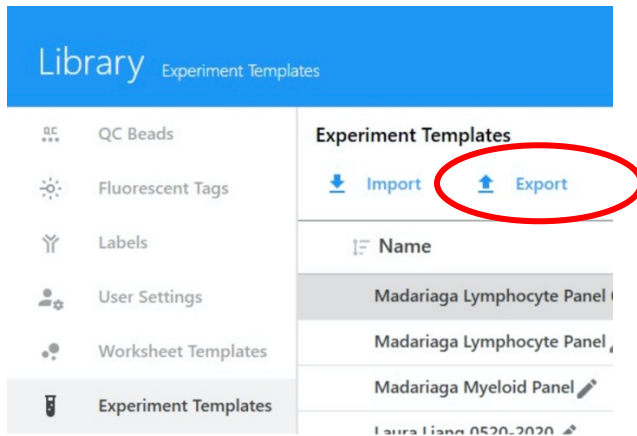

3. Save your template to the server or your preferred cloud storage. Saving will create a zip folder.

To import your saved template

1. In the acquisition tab, select import

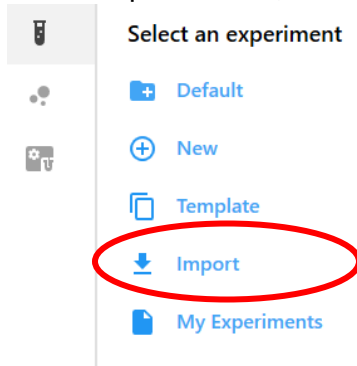

2. Locate your template (the zip folder) on the computer
3. A new experiment will automatically open from the experiment and you can run your samples
 - a. The template will now be saved in the template window
 - b. Don't forget to rename the new experiment to include your name and date!