

# IMPORTANT! Chamber Slide Users BEWARE!

You can *NOT* image the far wells for some designs on confocal systems with universal holders.

The objective will HIT THE HOLDER TABS if you try to focus, either breaking the lens and/or the stage.


**YOU WILL HAVE TO PAY FOR REPAIRS**

(\$1,200, \$6,000, or \$12,000)




Here's the problem. RED ARROWS show objective hits the support preventing FOCUS in yellow-tagged wells

OK wells in green, dangerous ones in red!

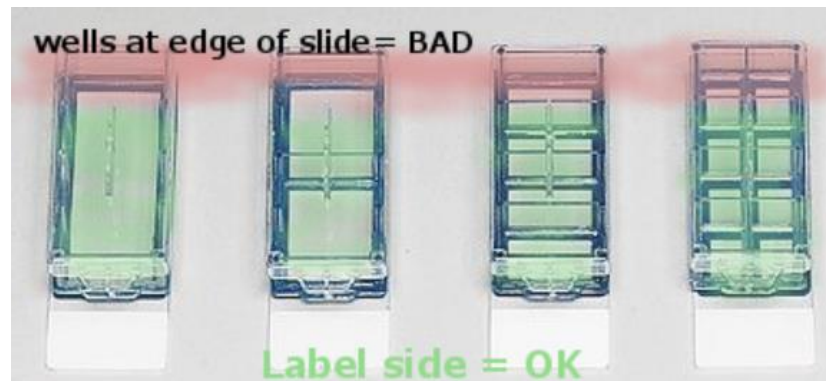

Note end tabs-- all are OK!

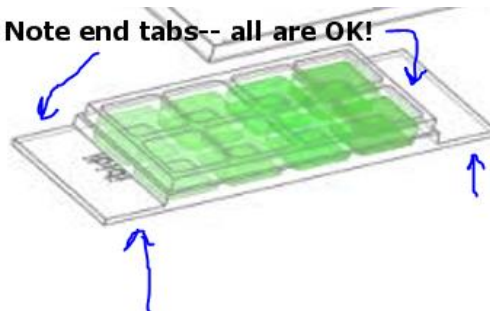

Coverslip slides:

